


FEMTOCONTROL

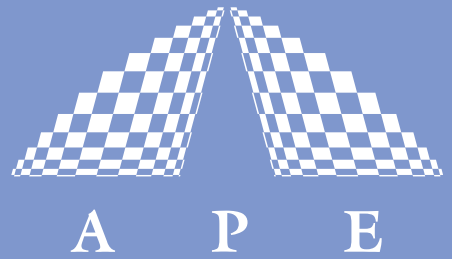

PULSE COMPRESSOR / PRE-CHIRPER



FemtoControl is a compact motorized pulse compressor for optimization of the duration of femto-second laser pulses in the spectral range of Ti:Sapphire lasers*.

Femtosecond pulses which pass through any kind of optical material experience dispersion that can lead to immense temporal broadening of the ultrashort pulses. Thus, the experimental conditions are changed and degraded.

FemtoControl compensates material dispersion by applying the inverse dispersion (or chirp) to the pulse. This inverse dispersion is generated by a pair of prisms on motorized translation stages allowing continuous adjustment of the pulse length.


Ideal for Multi-Photon-Microscopy

Wide range of dispersion

User-friendly adjustment and operation

Zero - GVD function

Automatic prism drive corresponding to stored GVD-curves

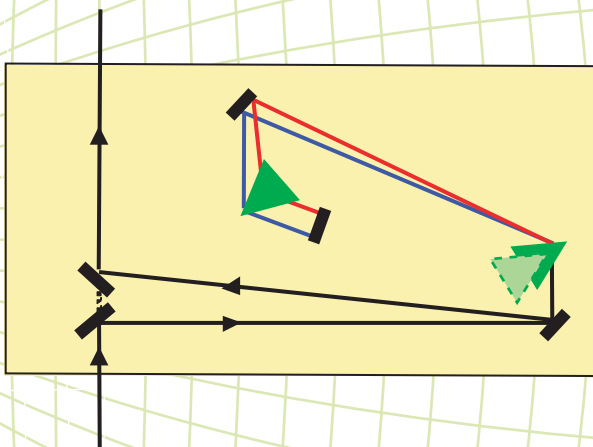

*For other laser systems contact APE GmbH

Ultrafast Pulse Diagnostics

Wavelength Conversion

Pulse Management

Acoustooptics

Your Partner in Ultrafast

FEMTOCONTROL


SPECIFICATIONS

Wavelength ranges	680 ... 1080 nm
Max. beam diameter	4 mm
Input polarization	Horizontal (polarization rotator optional)
Minimum laser repetition rate	Any
Max. bandwidth (for complete compensation)	12.5 nm @ 800 nm (\approx 75 fs transform limited Gaussian pulse; higher bandwidth on request)
Transmission	>90% @ 800 nm
Dispersion range	min max
	700 nm: 0 ... - 23000 fs ²
	800 nm: 0 ... - 13000 fs ²
	900 nm: 0 ... - 8000 fs ²
	(more dispersion on request)
Additional internal beam path	1.67 m

CONTROL OPTIONS

Remote control unit or control via autocorrelator **Carpe**

DIMENSIONS (in mm)

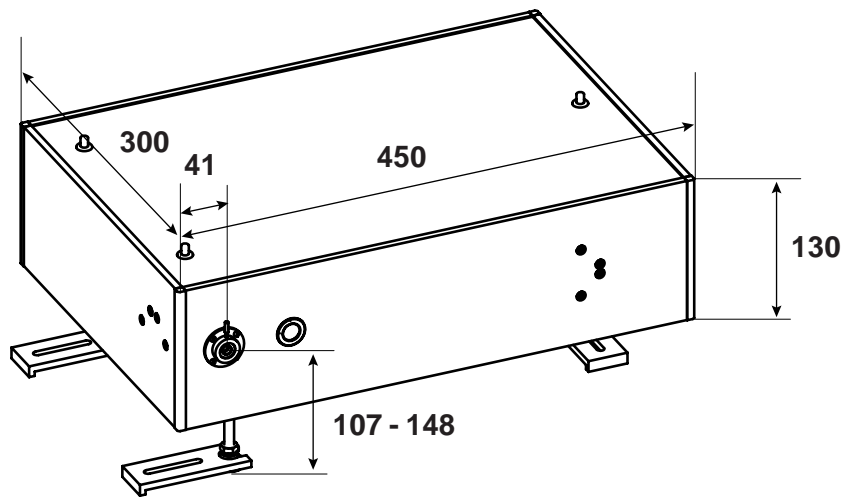

Application Example:

A transform limited Gaussian input pulse of 100 fs at 800 nm sent into a multi-photon microscope experiences a dispersion of 13000 fs² from the glass of the microscope lens system. That broadens the pulse to 380 fs at the sample position. **FemtoControl** in front of the microscope is capable of compensation back to 100 fs pulses at the sample.

At 700 nm a broadening of up to 645 fs, and at 900 nm of up to 245 fs can be compensated, respectively, for a 100 fs transform limited pulse.

Distributors

see APE website www.ape-berlin.com

APE GmbH Plauener Straße 163-165 Haus N / 13053 Berlin Germany
Phone +49.30.986.01130 Fax +49.30.986.011333 / Web www.ape-berlin.com Email ape@ape-berlin.de

APE follows a policy of continued product improvement. Therefore, specifications are subject to change without notice.